



Product: Anatomical So-Soft Fluid

Purpose: To preserve the soft tissues of the donor while maintaining a life-like texture and feel.

Useful length of time: 10 days without refrigeration. 6 weeks with refrigeration

How to use:

Mix ASSF by adding DI Water to a 2 water : 1 ASSF solution. The less water used the firmer the donor tissues will become. First time users follow this instruction before adjusting water volume on your first donor. You may add dye if you like.

Arterial Inject ASSF as you would a typical anatomical donor. Do not over distend the donor. Be sure that the vascular system runs clear before closing/tying off artery.

Allow donor to set in refrigeration for 24 hours prior to putting into lab.

Re-Order by emailing customerservice@trinityfluids.com