

# TRINITY FLUIDS, LLC

All Trinity Fluids, LLC brands are designed to promote positive results, if used properly. You can choose to not deviate from our instructions or you can experiment with it based on your level of experience and desired personal look and feel you desire.

## Usage instructions for Cadaver Hydrating Fluid Eco Blend, CHF-WS

- Mix 1 **CHF-WS** into 4 H<sub>2</sub>O. Blend continually for several minutes to ensure complete homogenous solution.
- Pour into spray.
- When ready to put cadaver away for extended periods apply topically **CHB-WS** all over areas in which you would like to maintain a soft and pliable condition. Do not over do it.
- Pour a half gallon of mixed **CHF-WS** into a bucket. Use cheese cloth or non cotton cloth to soak in solution. Twist and Squeeze out so cloth is not dripping with solution. Lay the cloth over the cadaver extremities and torso to increase hydration.

### **Optional additives to solution.**

- To KILL bacteria & excessive mold growth **add 6 ounce of SANISOL 7** to every one gallon of mixed solution.